

---

## Protein Block solution

#Cat: NB-23-00179-1

Size: 100ml (ready-to-use)

#Cat: NB-23-00179-2

Size: 18ml (ready-to-use)

### Intended Use

Protein Block Solution is a universal blocking reagent for reducing non-specific background staining in immunolabeling procedures. Its blocking capacity is superior to conventional blocking method. It eliminates the handling of animal serums in the laboratories. The solution is proven effective for immunohistochemistry, western blotting and ELISA procedures. It can be used as a diluting buffer for primary antibodies.

### Components

---

Contains 0.1%  $\text{NaN}_3$

### Storage

---

Store at 2-8°C.

### Recommended Protocol

---

1. Prior to application of primary antibody apply Protein Block Solution on tissue slides and incubate for 10 minutes at room temperature
2. Blot off Protein Block Solution (do not rinse) and add primary antibody

### Precaution

---

Wear gloves and take other necessary laboratory safety procedure.

**For Research Use Only. Not for Diagnostic or Therapeutic Use.**