# RT-MIS

ENGLISH .....	2
FRANCAIS.....	23
DEUTSCH.....	44
ESPAÑOL .....	65
ITALIANO .....	86

ENGLISH

## GENEXPATH RT-MIS software user guide

### User precautions.

*In vitro* diagnostic medical device according to Directive (EU) 98/79/EC

For *in vitro* diagnostic use

It is for professional use only.

Read all information in this user guide before use.

**Distributed by:**

**CliniSciences Group**

This document and its content are the property of GENEXPATH. They are only intended for contractual use by customers for the product described herein, excluding any other use. This document and its content may not be used or distributed for any other purpose, nor shared, disclosed or reproduced in any way without the prior written permission of GENEXPATH. GENEXPATH does not grant any license through this document.

The instructions given in this document must be strictly and explicitly followed by qualified and trained personnel to ensure appropriate use of the product described herein. All content in this document must be thoroughly read and understood before using the described product.

© 2024 Genexpath. All rights reserved.

ENGLISH.....	2
Foreword - please read carefully.....	5
About this guide.....	5
Product description .....	5
General information .....	5
GENEXPATH LymphoSign .....	5
GENEXPATH SarcomaFusion .....	5
GENEXPATH LymphoTranscript .....	6
System requirements.....	6
Home page.....	7
Analyses .....	8
Start a new analysis .....	8
Results.....	10
Processing details .....	10
LymphoSign Test .....	<b>Erreur ! Signet non défini.</b>
SarcomaFusion Test.....	<b>Erreur ! Signet non défini.</b>
LymphoTranscript Test .....	<b>Erreur ! Signet non défini.</b>
Specificities .....	19
Performances.....	19
Repeatability .....	19
LymphoSign Test .....	19
Random forest.....	19
Analytical specificity .....	19
Analytical sensitivity.....	19
Graph representation.....	19
SarcomaFusion Test .....	20
Associated Bibliography .....	21
Filtering non-specific fusion .....	20
Display of the fusions .....	20
SarcomaFusion and LymphoTranscript Tests .....	21
Internal control.....	21
Correction of hopping index.....	22
Notes.....	22
Reference.....	22

## Foreword - please read carefully

**GENEXPATH RT-MIS is a comprehensive bioinformatics solution which includes different data processing algorithms. It should be used critically by professionals in human genetics. Although GENEXPATH strives to offer high-quality software, it cannot guarantee the accuracy of the information and predictions it provides.**

## About this guide

This guide describes how to use the software GENEXPATH RT-MIS, Version 04, published in August 2024.

## Product description

### General information

After completing the *in vitro* protocol of a Genexpath test and sequencing, the software **GENEXPATH RT-MIS** allows the loading of sequencer files in FASTQ format, as well as the list of barcodes used during testing.

This software carries out demultiplexing to assign sequences to each sample. It then precisely identifies the gene expression markers and quantifies them.

It assesses the sequencing quality of each sample by quantifying the number of reads identified and the number of UMI (unique molecular identifiers) detected.

For each sample, it generates a report to help the user analyse the data. This data can be downloaded.

### GENEXPATH LymphoSign

Through the use of LymphoSign tests, the software helps analyse results obtained after completing the *in vitro* protocol and sequencing. The analysis is based both on quantification of gene marker expression, but also the presence or absence of mutations and chromosome translocations.

The artificial intelligence classification model was trained with a database containing over 3000 analyses.

If this database were to change, users will be kept informed via the RT-MIS platform.

The expression level of each marker is normalised in relation to the average gene expression of all markers studied, allowing a comparison of analyses.

### GENEXPATH SarcomaFusion

Through the use of SarcomaFusion tests, the software helps analyse results obtained after completing the *in vitro* protocol and sequencing. RT-MIS provides information on the presence or absence of chromosome translocations amongst 140 determined fusions associated with

sarcomas. It also establishes a link with bibliography references associated with detected fusions.

The bibliography used was created from Mitelman databases [1] and PubMed research for all fusion transcripts associated with sarcomas. It is non-exhaustive and specific to sarcomas.

### **GENEXPATH LymphoTranscript**

Through the use of LymphoTranscript tests, the software helps analyse results obtained after completing the *in vitro* protocol and sequencing. RT-MIS provides information on the presence or absence of chromosome translocations by studying 69 genes associated with T-cell non-Hodgkin's lymphomas.

### **GENEXPATH CustomPanel**

In the context of the use of CustomPanel assays, the software allows the analysis of the results obtained after the *in vitro* protocol and sequencing have been performed. The raw data and expression level of each marker are provided. The expression level of each marker is normalized in relation to the average gene expression of all the markers studied, thus allowing the comparison of the analyses with each other.

### **System requirements**

The software RT-MIS is directly accessible through the user space at <https://connect.genexpath.com>. Access is restricted to user accounts authorised by Genexpath.

To get an account, send an email to the following address: [support@genexpath.com](mailto:support@genexpath.com).

An internet connection is required. The software is compatible with the browsers Mozilla Firefox, Microsoft Edge and Google Chrome. It does not require the installation of additional tools.

## Home page

After logging in to the user space, RT-MIS is accessed by clicking on the RT-MIS button on the left-hand bar, then on “RT-MIS” (Figure 1):

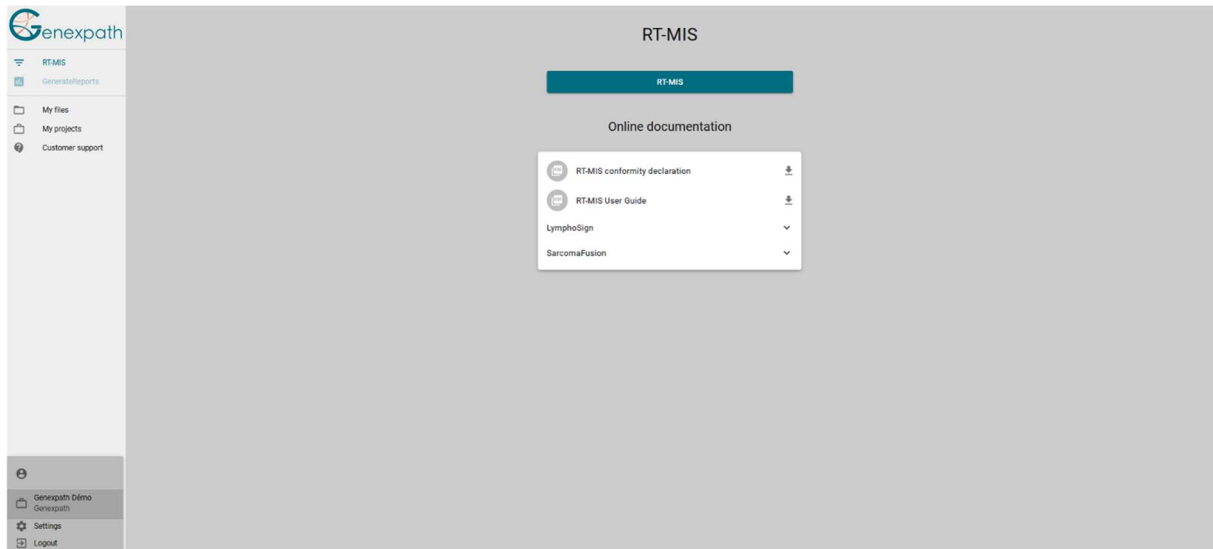
Figure 1: User space page

The home page is shown (Figure 2):

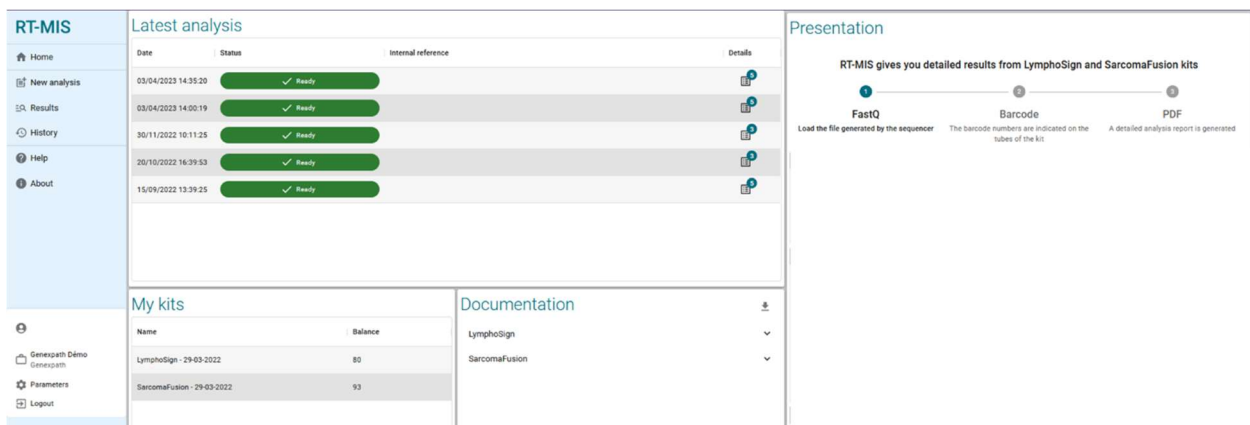
Figure 2: RT-MIS home page

It contains several windows:

- **“Latest analysis”** with the latest analyses completed, their status and access to analysis detail by clicking on the corresponding icon . The number indicated in the top right corresponds to the number of samples processed (and therefore deducted from the corresponding kit balance).
- **“Presentation”** briefly summarises how RT-MIS works, as well the latest platform updates/news.

- **“My kits”** contains information on the kits held by the centre to which the user is linked, and the remaining analysis balance.
- **“Documentation”** contains a link to Genexpath product user guides.
- **“Updates”** contains last information of updates.

On the left, the main menu (top) allows you to access:

- On the home page
- New analysis launch
- Display of results
- kits ‘history
- Interactive help (for the home page and the “new analysis” page)
- About

The main menu (bottom) indicates the login with which you are connected and allows:

- To return to the user area
- To choose to receive an email at the end of an analysis
- Sign out.

## Analyses

### Start a new analysis

- 1) In the menu on the left, click on **“New analysis”** (Figure 3).

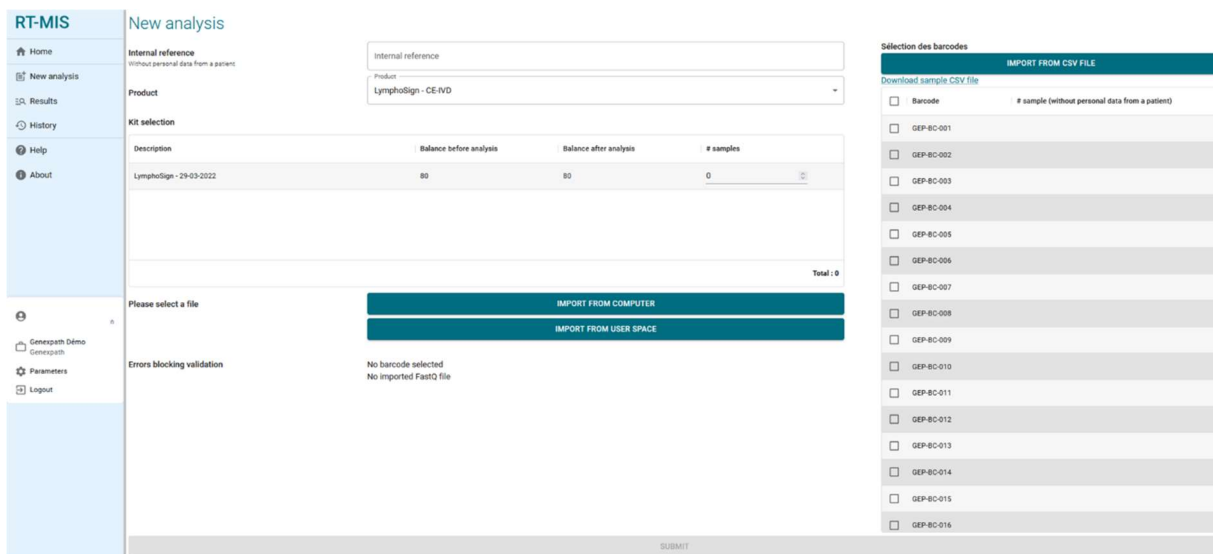
Figure 3: “New analysis” page

- 2) Indicate an internal reference to the analysis with any patient data information (figure 4)
- 3) Select in the dropdown menu “product”, the used product.
- 4) In “kit selection”, indicate the number of analyses that you need to count down per kit (according to the batch number used in the “*in vitro*” part).
- 5) Select a FastQ file to analyse by clicking on:
  - **“IMPORT FROM COMPUTER”** for a file saved on the computer, or
  - **“IMPORT FROM USER SPACE”** for a file already imported to the user space.

The FASTQ file must respect the naming convention of Illumina sequencers.

MiSeq example: X\_S1\_L001\_R1\_001.fastq.gz

The software checks the format and integrity of the FastQ file(s) before proceeding to the analysis. This check lasts from a few seconds to a few minutes, depending on the file size.

- 6) Tick (right side of the screen) the barcodes used for the analysis (numbers printed on the tubes).

For each selected barcode, an analysis will be deducted.

You can add an identifier for each selected barcode.

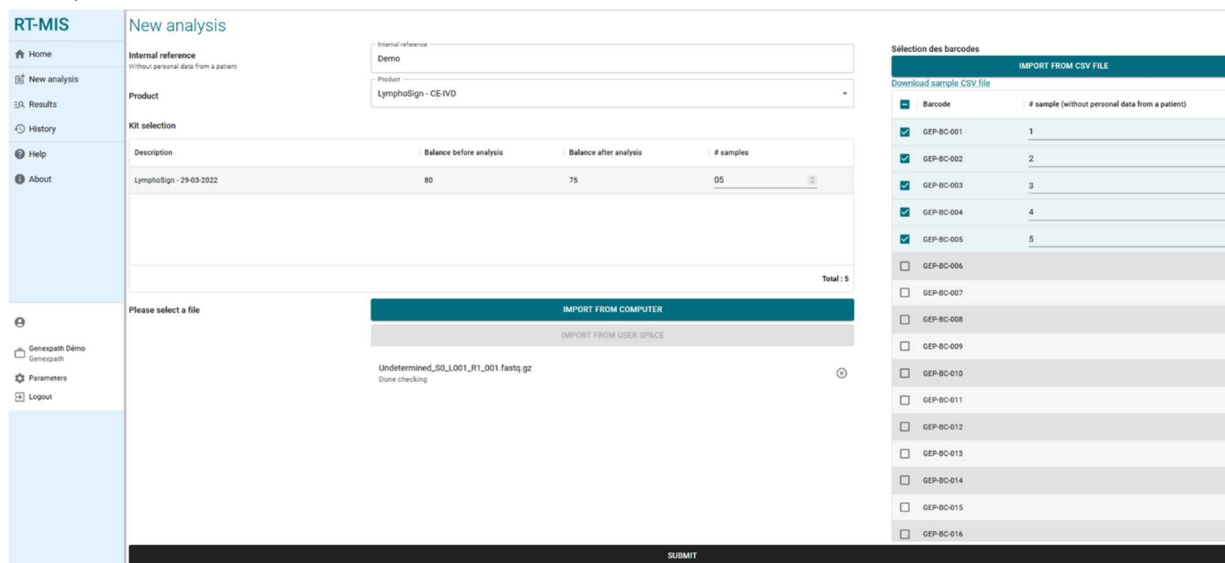
No patient data can be entered.

You can also have the option of downloading directly a CSV file with its information from the CSV file example.

Issues that could block the validation are:

- Absence of FASTQ file
- Barcodes selection mismatch

- 7) Click on **“VALIDATE”**.

Description	Balance before analysis	Balance after analysis	# samples
LymphoSign - 29-03-2022	80	75	05

Barcode	# sample (without personal data from a patient)
<input checked="" type="checkbox"/> GEP-BC-001	1
<input checked="" type="checkbox"/> GEP-BC-002	2
<input checked="" type="checkbox"/> GEP-BC-003	3
<input checked="" type="checkbox"/> GEP-BC-004	4
<input checked="" type="checkbox"/> GEP-BC-005	5
<input type="checkbox"/> GEP-BC-006	
<input type="checkbox"/> GEP-BC-007	
<input type="checkbox"/> GEP-BC-008	
<input type="checkbox"/> GEP-BC-009	
<input type="checkbox"/> GEP-BC-010	
<input type="checkbox"/> GEP-BC-011	
<input type="checkbox"/> GEP-BC-012	
<input type="checkbox"/> GEP-BC-013	
<input type="checkbox"/> GEP-BC-014	
<input type="checkbox"/> GEP-BC-015	
<input type="checkbox"/> GEP-BC-016	

Figure 4: Information on the user analysis

RT-MIS opens the **“Results”** page.

## Results

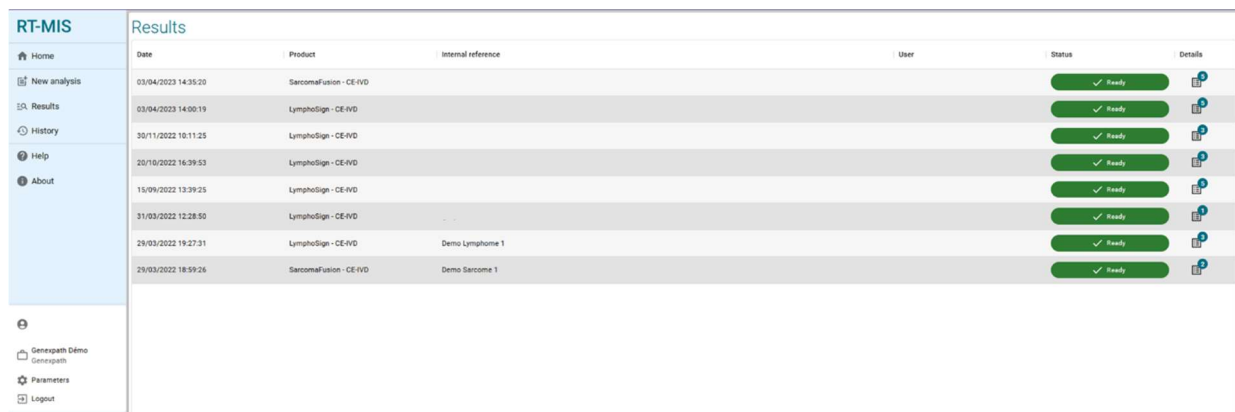
On the left side menu, click on item **“Results”** (figure 5).

This page offers access to new and previous results.

RT-MIS indicates the analysis status: **“pending processing”**, **“processing”** or **“processing complete”**.

The results will be available once processing is complete. Processing can last several minutes.

To access analysis details, click on the corresponding icon .

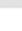
Date	Product	Internal reference	User	Status	Details
03/04/2023 14:35:20	SarcomaFusion - CE-IVD			✓ Ready	
03/04/2023 14:00:19	LymphoSign - CE-IVD			✓ Ready	
20/11/2022 10:11:25	LymphoSign - CE-IVD			✓ Ready	
20/10/2022 16:39:53	LymphoSign - CE-IVD			✓ Ready	
15/09/2022 13:39:25	LymphoSign - CE-IVD			✓ Ready	
31/03/2022 12:28:50	LymphoSign - CE-IVD			✓ Ready	
28/03/2022 19:27:31	LymphoSign - CE-IVD	Demo Lymphome 1		✓ Ready	
29/03/2022 18:59:26	SarcomaFusion - CE-IVD	Demo Sarcome 1		✓ Ready	

Figure 5: “Results” page

It is possible to switch pages in the bottom right of the window, depending on the number of analyses completed.

## Processing details

### All tests

This page offers a summary of the analysis characteristics (Figure 6).

- The status
- The centre
- The project
- The internal reference
- The product used
- The user
- The FASTQ file save date
- The FASTQ file processing date
- The kit used and the number of analyses used on it.

Process detail

Status	✓ Ready	
Center	Genexpath	
Laboratory	Genexpath Rouen	
Internal reference		
Product	LymphoSign - RUO	
User		
Record date	05/08/2024 13:19:39	
Process date	05/08/2024 13:20:02	
Kit	Kit	Analysis consumed on the kit
	LymphoSign	11

Figure 6: "Processing details" page

Quality metrics are represented for the entire treatment (Figure 7):

- Genexpath barcodes: Percentage of reads associated with Genexpath libraries on the entire FASTQ file loaded.
- Percentages of barcoded reads according to its barcode
- Representation of the reads of the loaded FASTQ file

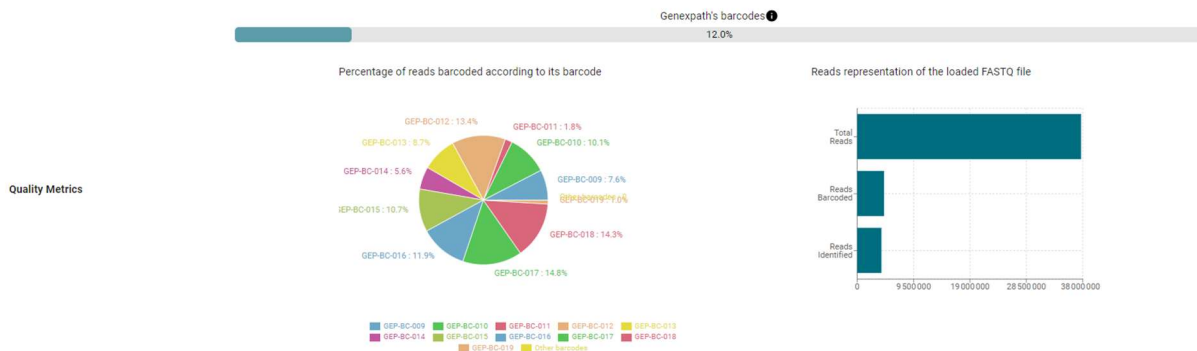
Figure 7: Quality metrics

For each sample tested, you will find the results of your analysis with a reminder of the barcode used and the sample reference, if applicable. This part allows you to observe the informativity for all the samples in the analysis (Figure 8).

Barcode	Sample	Informativity LymphoSign - RUO	Reports	Count data
<input checked="" type="checkbox"/> GEP-BC-009		88.5%		
<input checked="" type="checkbox"/> GEP-BC-010		91.8%		
<input checked="" type="checkbox"/> GEP-BC-011		63.6%		
<input checked="" type="checkbox"/> GEP-BC-012		92.7%		
<input checked="" type="checkbox"/> GEP-BC-013		90.6%		
<input checked="" type="checkbox"/> GEP-BC-014		89.9%		
<input checked="" type="checkbox"/> GEP-BC-015		92.6%		
<input checked="" type="checkbox"/> GEP-BC-016		93.2%		
<input checked="" type="checkbox"/> GEP-BC-017		89%		
<input checked="" type="checkbox"/> GEP-BC-018		93.5%		
<input checked="" type="checkbox"/> GEP-BC-019		4.5%		

Download DOWNLOAD ALL DOCUMENTS DOWNLOAD ALL COUNT DATA

Figure 8: Access to analyses by barcode

## LymphoSign Test

It is possible to generate a customizable analysis report by clicking on .  
A report generation page opens, by default on tab "results":

### Identification

Sample		Further information	
Barcode	GEP-BC-013	User	
# sample		Process date	05/08/2024 13:20:02
Center	Genexpath	Application version	
Laboratory	Genexpath Rouen	Internal reference	
Kits	LymphoSign		

### Quality control

Number of reads	Number of UMI	Informativity in %	Amplification factor
396068	33905	90.6	11.7

### Expression signature

Predictions must be correlated with histopathological analysis.  
Barcode: GEP-BC-013  
Predicted GEP class: **DLBCL ABC**  
Additional Features: **MYD88L265P**  
Quality score: **33905** (≥ 5000 - pass)

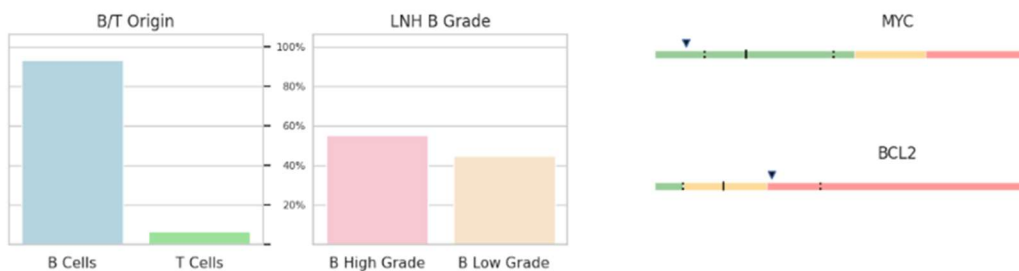
Figure 9: Example report for LymphoSign test

By default, the report contains the following information:

- Analysis identification: reminder of analysis characteristics
- Quality control: number of reads and UMI (Unique Molecular Identifiers), the informativity (number of junction L probes – R waited probes / total number of reads \*100) and the amplification factor.
- The expression signature: barcode, predicted class, optional additional feature, quality score, as well as a set of graphs detailing the inferences made to predict the class of non-Hodgkin's lymphoma.

Name	Description
<b>Barcode</b>	Reminder of barcode used for the analysis
<b>Predicted GEP class</b>	Class based on AI-predicted gene expression profile
<b>Additional Features</b>	Additional information
<b>Quality score</b>	Quality score based on number of UMI

You can also choose which items you will eventually see appear in the report using the "Add Item" buttons depending on how many charts you will be able to add. In addition, you can view the report before downloading by clicking on "view PDF". The graph is downloaded by clicking on "download PDF".

The additional elements are:

- A customizable graph (Figure 10) representing the level of expression of several genetic markers of interest according to the position of your sample (red dot) as well as:
  - o the entire database
  - o One of the two classes LNH\_B or LNH\_T
  - o Of a sub-entity

Custom graph

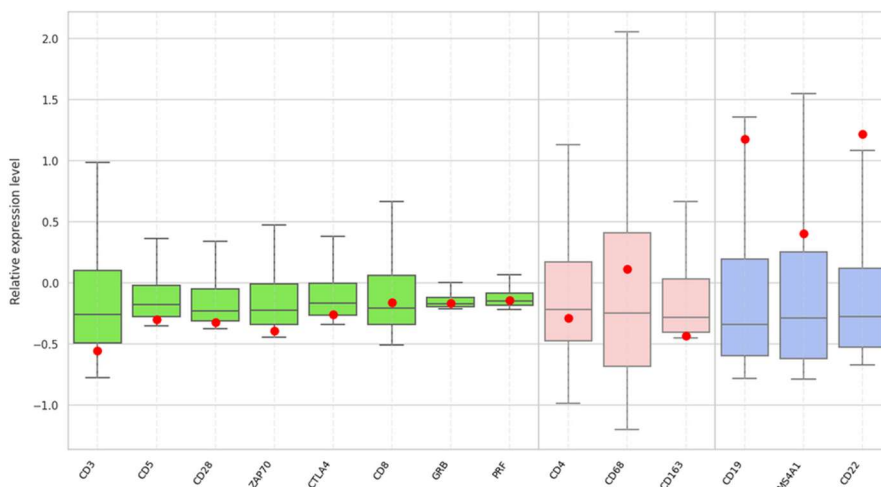
Figure 10: Generation of custom graph

You can at the top right (Figure 11):

- o Choose the number of classes wished by clicking on “add a class”
- o Determine the name of each class
- o Choose the markers of interest by clicking on the pencil of the corresponding class

Edit class

Class name

---

CD3 CD5 CD28 ZAP70

CTLA4 CD8 GRB PRF

Associated color

SUBMIT

*Figure 11: Determination of markers for a given class*

- Predefined graphical representations with the comparison of the level of expression (Figure 12). The reference expression levels are displayed according to the chosen class and the sample analyzed is represented by a black diamond or a red dot depending on the tab. It is possible to compare gene expression levels.

Tab	Description
<b>Comparison</b>	Comparison of expression profiles between two LymphoSign signature classes (the predicted one and another chosen by the user)
<b>B vs T</b>	Specific B and T lymphocyte markers
<b>IG</b>	Different immunoglobulins
<b>MUTS</b>	Specific markers such as point mutations
<b>DLBCL</b>	Specific markers of diffuse large B cell lymphomas (DLBCL)
<b>PTCL</b>	Specific markers of peripheral T-cell lymphomas
<b>Other</b>	Specific markers of different lymphomas
<b>Prognosis</b>	Markers which could affect patient prognosis

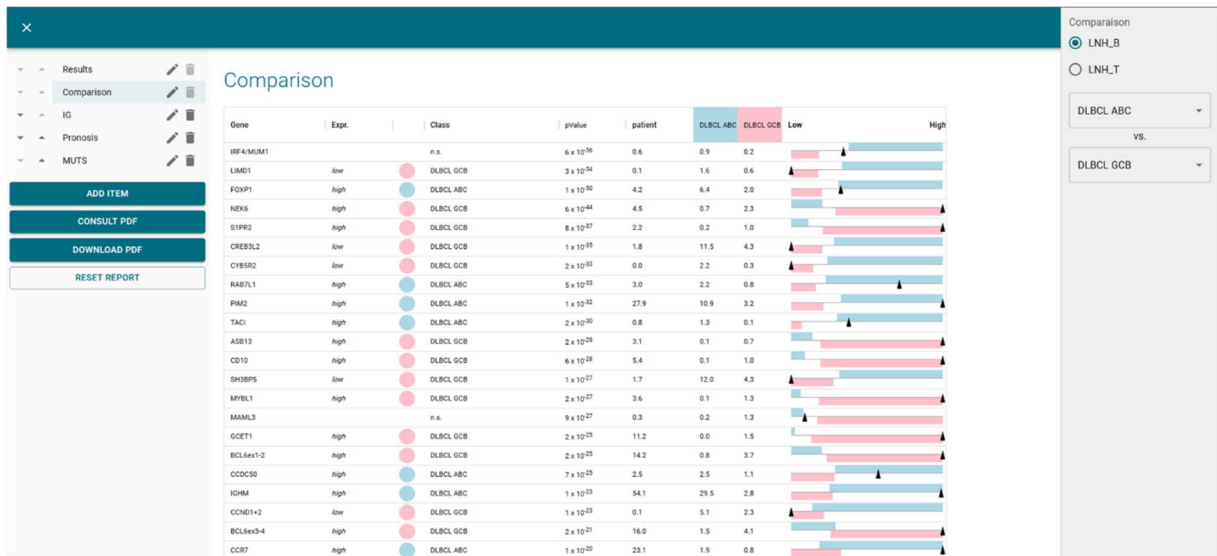
Figure 12: Example of advanced interpretation with comparison between PMBL and GCB

Only the first 30 genes of the two-class comparison are displayed.

Each of these comparisons can be personalised using the right-hand menu, through which you can select the type of classes to compare (Figure 13):

- The entire database
- one of the two classes LNH\_B or LNH\_T
- A sub-entity

Only the first 30 genes of the comparison of 2 classes will be display.

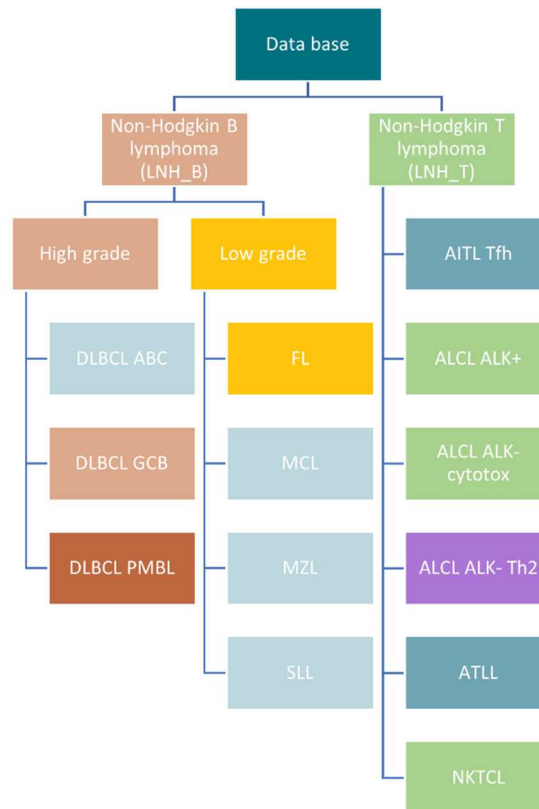
Figure 13: Representation of different classes predicted by RT-MIS

You may also download in “**management details**” page:

- For each sample: the CVS file of all raw and relative data of by clicking on 
- The CSV file with all raw and relative data by clicking on “**DOWNLOAD ALL COUNTING DATA**”
- All pdf report or CSV files can be downloaded by clicking on “**ALL REPORT**”

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By clicking on  icon of barcode of interest, you can also generate a report (figure 14) with or without the associated bibliography, consult and download it.

Identification

Sample		Further information	
Barcode	GEP-BC-002	User	
# sample		Process date	01/08/2024 13:37:01
Center	Genexpath	Application version	
Laboratory	Genexpath Rouen	Internal reference	
Kits	SarcomaFusion		

Quality control

Quality control SarcomaFusion		Quality control GAPDH	
Number of reads	189638	Number of reads	187109
Number of UMI	2954	Number of UMI	11767
Amplification factor	64.2	Amplification factor	15.9
Informativity	85.4%	Informativity GAPDH	95.7%

Results

Fusion	Number of reads	Number of UMI	Amplification factor	References
✓ PAX3_7E7G-FOXO2D	158329	1163	136.1	<a href="#">7 references</a>

Fusion : PAX3\_7E7G-FOXO2D

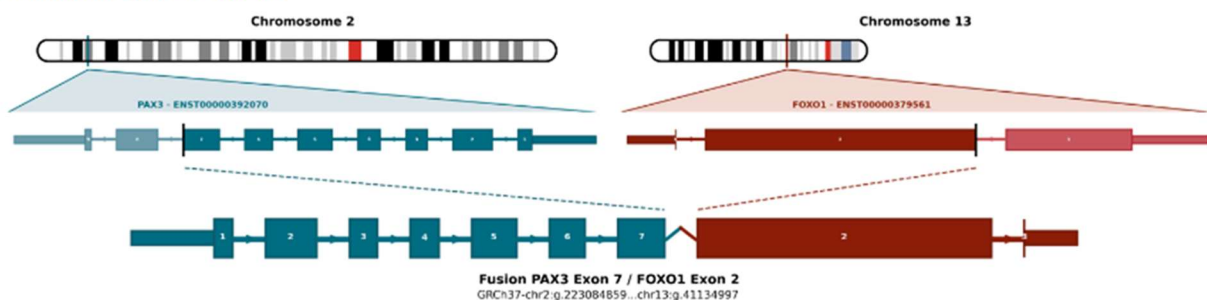
Figure 14: "Report" page for the SarcomaFusion test

Information available:


- Result with
  - Summary of the analysis characteristics
  - Quality control
  - number of reads and UMI for the sample of interest.
  - The informativity in % as well as the amplification factor.
  - number of reads and UMI for the GAPDH internal control
  - The informativity in % as well as the amplification factor for internal control

- The results of the fusion may have been recognized
- the bibliography references if the fusion is associated to a sarcoma or a T-cell lymphoma.
- The raw data obtained by RT-MIS, without filtration of the number of UMI, with a graphic representation of the detected fusion depending on the number of reads or UMI and a summarize table.
- A graphic representation of the raw data

You may also download in “**management details**” page:

- For each sample: the CVS file of all raw and relative data of by clicking on 
- The CSV file with all raw and relative data by clicking on “**DOWNLOAD ALL COUNT DATA**”
- All pdf report or CSV files can be downloaded by clicking on “**DOWNLOAD ALL REPORT**”

Fusion transcript notation in RT-MIS is as follows (Figure 15):

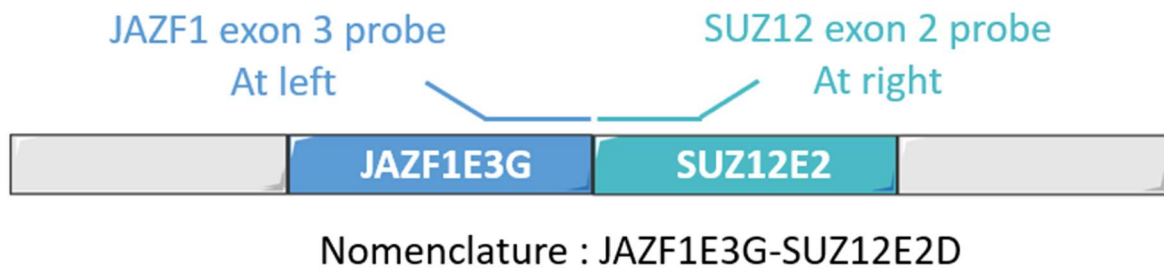*Figure 15: Schematic representation of naming system for fusion transcripts*

By clicking on “n references”, the bibliography details open in a new page (Figure 16).

#### Bibliographic report

Name	References
JAZF1/SUZ12	<a href="#">Novel EPC1 gene fusions in endometrial stromal sarcoma.</a> Genes Chromosomes Cancer
JAZF1/SUZ12	<a href="#">Cytogenetic and molecular genetic analyses of endometrial stromal sarcoma: nonrandom involvement of chromosome arms 6p and 7p and confirmation of JAZF1/JJAZ1 gene fusion in t(7;17).</a> Cancer Genet Cytogenet
JAZF1/SUZ12	<a href="#">Molecular analysis of the JAZF1-JJAZ1 gene fusion by RT-PCR and fluorescence in situ hybridization in endometrial stromal neoplasms.</a> Am J Surg Pathol
JAZF1/SUZ12	<a href="#">Extrauterine endometrial stromal sarcoma with JAZF1/JJAZ1 fusion confirmed by RT-PCR and interphase FISH presenting as an inguinal tumor.</a> Virchows Arch
JAZF1/SUZ12	<a href="#">A neoplastic gene fusion mimics trans-splicing of RNAs in normal human cells.</a> Science
JAZF1/SUZ12	<a href="#">Frequent fusion of the JAZF1 and JJAZ1 genes in endometrial stromal tumors.</a> Proc Natl Acad Sci U S A

*Figure 16: “Bibliography references” page*

## Specificities

### Performances

#### *Repeatability*

The repeatability of the RT-MIS software is defined as its ability to obtain the same results. The same FASTQ analyzed in triplicate with 15 samples for each of two tests was studied. The RT-MIS software has a repeatability of 100%. The same FASTQ file containing several samples gives:

- for a LymphoSign test: the same classification as well as the same expression levels for each of the markers
- for a SarcomaFusion test: the same result (presence or absence of a fusion) as well as the associated bibliography.

#### **LymphoSign Test**

#### *Random forest*

Normalization of expression data is a step aimed at making samples sequenced under different experimental conditions comparable (quantity and quality of RNA, type of sequencer, number of reads generated, etc.). In order to erase these experimental biases, the raw counting data is divided by the median of the counts of a fixed list of markers.

From the normalized expression data, RT-MIS classifies the sample by Random Forest (RF). This classification is based on the application to the expression data of 10 RF classification models from the Machine Learning software from the database of nearly 3000 cases. The different probabilities of these classifiers make it possible to identify different signatures and classes described in Figure 13.

#### *Analytical specificity*

A study was carried out on a larger scale with 283 cases (Bobee et al., Blood Cancer Journal, 2020). The objective of this study was to demonstrate the relevance/power of the Random Forest algorithm allowing classification by comparing the results with another classification technique. 94.3% of the results were concordant. Thanks to the algorithm, some of the cases were later reclassified.

#### *Analytical sensitivity*

The artificial intelligence was trained on a database of nearly 3,000 cases. The prediction is optimal when there are at least 5000 UMI.

#### *Graph representation*

The level of expression of each of the markers corresponds to the raw count data divided by the median of the counts of a fixed list of markers.

The data is represented in the form of boxplots. They correspond for a given marker to the reference values of the database according to the entire database, of one of the two classes LNH\_B or LNH\_T or of a sub-entity.

Each boxplot is organized as follows:

- The central value of the graph is the median
- The edges of the rectangle are the quartiles (Q1 at the bottom and Q3 at the top)
- The endpoints of the boxplot are calculated using 1.5 times the interquartile space (the distance between the 1st and 3rd quartiles).

## SarcomaFusion Test

### *Filtering non-specific fusion*

The background noise is established by a list of non-specific fusions which is filtered to facilitate the understanding of the results. It was created from the results obtained from 5 tested samples whose expected fusions were known. Non-specific fusions were detected in at least 2 samples.

This list is as follows:

SS18E10G-SS18E3D	FUSE5G-STAT6E17D	RAD51BE3G-TFE3E7D
NAB2E3G-DDIT3E3D	HEY1E4G-CREB3L2E6D	FUSE9bG-CAMTA1E9D
PAX3_7E7G-TFE3E3D	KIF5BE24G-CAMTA1E9D	CTNNB1E1G-NTRK1E16D
ZC3H7BE12G-RELA11D	RBPMSE5G-EGFE17D	COL1A1E37G-STAT6E17D
TEAD1E8G-ETV1E4D	EWSR1E15G-ALKE20delD	CTNNB1E1G-NTRK3E18D
CITED2E2G-ERGE10D	LIFRE1G-NFATC2E3D	
HEY1E4G-NTRK1E10D	TFGE4G-NTRK3E11D	

These non-specific fusions do not appear in the PDF report, nor in the CSV file which can be downloaded.

### *Display of the fusions*

The pairs of probes detected correspond to potential genetic anomalies. They are compared to the bibliographic database. The objective is to determine whether the fusion detected has already been described in the literature in the context of sarcomas or not. Depending on this referencing, the merge is retained or filtered according to the algorithm below (Figure 17):

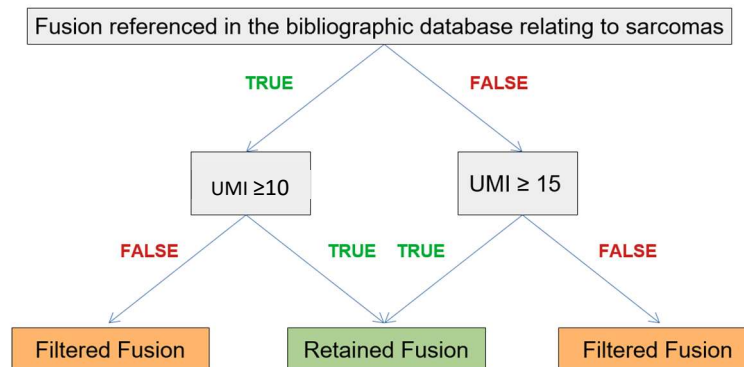
Figure 17: Algorithm of filtration representation

The retained fusion (or fusions) is returned as a result in the "Filtered data" part of the report and, if applicable, associated with its literature and the filtered fusions are entered in the "Raw data" part.

## LymphoTranscript Test

### *Filtration of non-specific fusions*

The background is established by a list of non-specific merges that is filtered to make the results easier to understand. It was created from the results obtained from 5 tested samples whose expected melts were known. Non-specific fusions were detected in at least 3 samples. This list is as follows: CD74E7G-DUSP22E2D, TPM3E8G-TPM3E3D.

These non-specific merges do not appear in the downloadable PDF report or CSV file.

### *Viewing Fusion transcripts*

The pairs of probes detected correspond to potential genetic abnormalities. They are compared with the bibliographic database. The objective is to determine whether the detected fusion has already been described in the literature in a T-cell lymphoma context or not. Based on this referencing, the fusion is kept above 10 UMI.

The retained fusion(s) is rendered as a result in the "Filtered data" section of the report and, if applicable, associated with its literature, and the filtered fusions are populated in the "Raw data" section.

## SarcomaFusion and LymphoTranscript Tests

### *Associated Bibliography*

The bibliographic database used was created from the Mitelman databases [1] and Pubmed searches for all the fusion transcripts associated with sarcomas. It is non-exhaustive and specific to sarcomas. It will be updated twice a year.

### *Internal control*

It is recommended to perform an internal control during the *in vitro* test using the same barcode and control barcode number for the same sample, as indicated in the SarcomaFusion and LymphoTranscript instructions for use. During computer analysis, RT-MIS combines the raw data from both barcodes to generate a single report.

The sample with the barcode must be sequenced as a normal sample with 100 000 reads. Nevertheless, its usage has no impact on the analysis countdown.

### *Correction of hopping index*

Hopping index is a phenomenon, during sequencing, of exchanges of barcodes leading to a wrong attribution of the sequence/sample couple. This phenomenon, in the case of the SarcomaFusion and LymphoTranscript tests, can lead to misinterpretation of the results.

To correct this technical effect, the UMIs assigned to each pair of left/right probes of the different sequenced samples are compared with each other. The UMI sequences shared by the same pair of probes between two barcodes of the same run are eliminated when they are not amplified.

## Notes

The company **GENEXPATH** does not permanently store results generated by the software **GENEXPATH RT-MIS**. The data must be downloaded directly after each analysis and saved by the user in their document management system.

However, analysis results are available for one year.

Information in this document is likely to change.

## Reference

[1] Mitelman Database of Chromosome Aberrations and Gene Fusions in Cancer (2021). Mitelman F, Johansson B and Mertens F (Eds.), <https://mitelmandatabase.isb-cgc.org>

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RT-MIS	
Nom	Référence
PAX3/FOXO1	Alveolar rhabdomyosarcoma of the head and neck region in older adults: genetic characterization and a review of the literature. Hum Pathol
PAX3/FOXO1	Cryptic insertion of 3'FOXO1 into inverted chromosome arm 2q in the presence of two normal chromosome 13s and 13 small interstitial duplications in a patient with alveolar rhabdomyosarcoma. Cancer Gene
PAX3/FOXO1	Molecular Cytogenetics Detect an Unbalanced t(2;13)(q26;q14) and PAX3-FOXO1 Fusion in Rhabdomyosarcoma With Mixed Embryonal/Alveolar Features. Pediatr Blood Cancer
PAX3/FOXO1	Molecular diagnostics in the management of rhabdomyosarcoma. Expert Rev Mol Diagn
PAX3/FOXO1	Clinicopathologic and Molecular Features of a Series of 41 Biphenotypic Sinonasal Sarcomas Expanding Their Molecular Spectrum. Am J Surg Pathol
PAX3/FOXO1	Early detection of the PAX3-FOXO1 fusion gene in circulating tumor-derived DNA in a case of alveolar rhabdomyosarcoma. Genes Chromosomes Cancer
PAX3/FOXO1	Alternate PAX3-FOXO1 oncogenic fusion in biphenotypic sinonasal sarcoma. Genes Chromosomes Cancer

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Process detail					
Status	<span style="color: green;">✓ Ready</span>				
Center	Genexpath				
Laboratory	Genexpath Rouen				
Internal reference					
Product	LymphoSign - RUO				
User					
Record date	05/08/2024 13:19:39				
Process date	05/08/2024 13:20:02				
Kit	<table border="1"> <thead> <tr> <th>Kit</th> <th>Analysis consumed on the kit</th> </tr> </thead> <tbody> <tr> <td>LymphoSign</td> <td>11</td> </tr> </tbody> </table>	Kit	Analysis consumed on the kit	LymphoSign	11
Kit	Analysis consumed on the kit				
LymphoSign	11				

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### Expression signature

Predictions must be correlated with histopathological analysis.

Barcode: GEP-BC-013  
 Predicted GEP class: **DLBCL ABC**  
 Additional Features: **MYD88L265P**  
 Quality score: **33905** (≥ 5000 - pass)

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### Identification

Sample	
Barcode	GEP-BC-002
# sample	
Center	Genexpath
Laboratory	Genexpath Rouen
Kits	SarcomaFusion

Further information	
User	
Process date	01/08/2024 13:37:01
Application version	
Internal reference	

### Quality control

Quality control SarcomaFusion	
Number of reads	189638
Number of UMI	2954
Amplification factor	64.2
Informativity	
85.4%	

Quality control GAPDH	
Number of reads	187109
Number of UMI	11767
Amplification factor	15.9
Informativity GAPDH	
95.7%	

### Results

Fusion	Number of reads	Number of UMI	Amplification factor	References
✓ PAX3_7E7G-FOXO2D	158329	1163	136.1	7 references

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
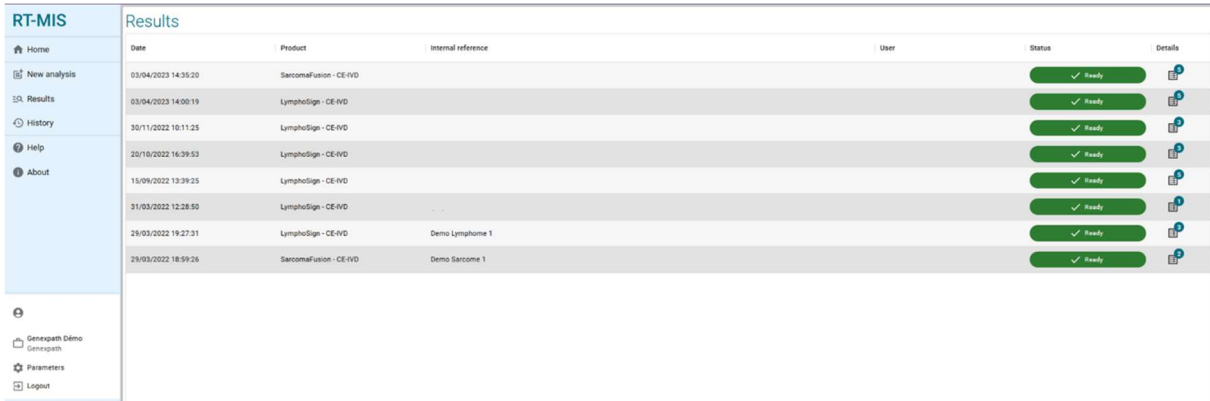
RT-MIS		Results				
Home	Date	Product	Internal reference	User	Status	Details
New analysis	03/04/2023 14:35:20	SarcomaFusion - CE-IVD			✓ Ready	
Results	03/04/2023 14:00:19	LymphoSign - CE-IVD			✓ Ready	
History	30/11/2022 10:11:25	LymphoSign - CE-IVD			✓ Ready	
Help	20/10/2022 16:39:53	LymphoSign - CE-IVD			✓ Ready	
About	15/09/2022 13:39:25	LymphoSign - CE-IVD			✓ Ready	
	31/03/2022 12:28:50	LymphoSign - CE-IVD			✓ Ready	
	29/03/2022 19:27:31	LymphoSign - CE-IVD	Demo Lymphome 1		✓ Ready	
	29/03/2022 18:39:26	SarcomaFusion - CE-IVD	Demo Sarcome 1		✓ Ready	

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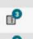Download  DOWNLOAD ALL DOCUMENTS  DOWNLOAD ALL COUNT DATA

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Company: CliniSciences AG
Company: CliniSciences Limited
Company: Biotrend Chemicals LLC